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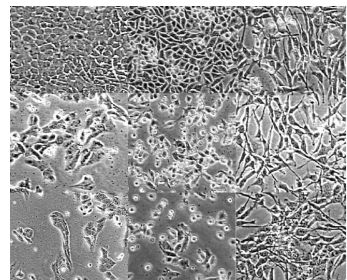
Dilution Calculator

Current concentration: 5.63×10^5 /mL Live cells

What is your desired cell concentration? x Live cells/mL Total cells

How many mL do you need?



Mix **1.56** mL of your cell solution with **10.44** mL buffer.

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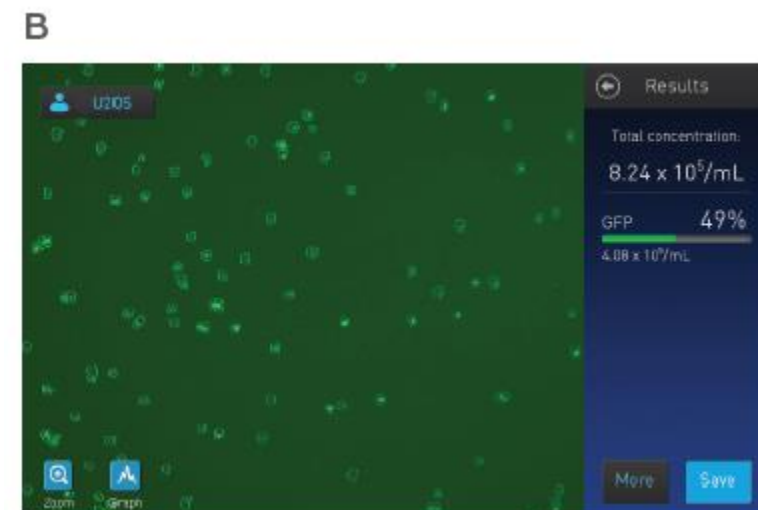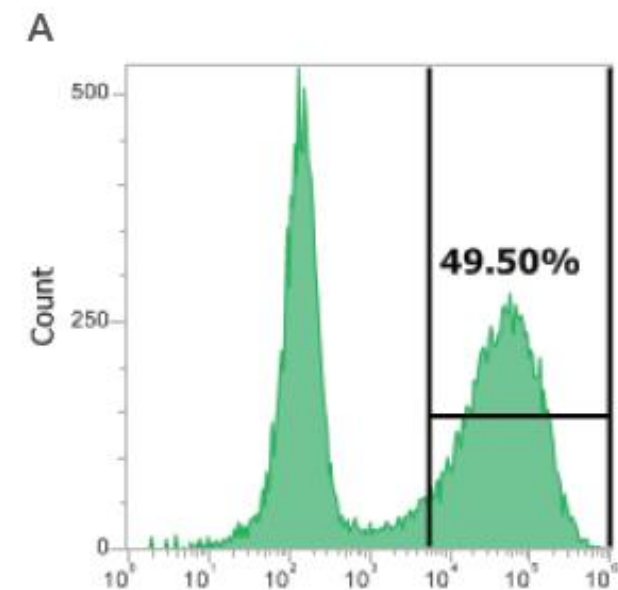For a complete list of available common and specialty light cubes, go to [thermofisher.com/evoslightcubes](https://www.thermofisher.com/evoslightcubes)

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invitrogen

Click the name of a light cube to view the list of compatible dyes

Commonly used light cubes

DAPI

TagBFP

CFP

GFP

YFP

RFP

Texas Red™

Cy5

EVOS Light Cube selection guide

Invitrogen™ EVOS™ Light Cubes integrate LEDs and high-performance filters in a novel illumination system that is designed to deliver precise control, minimal maintenance, and exceptional reliability. These LED cubes are compatible with the EVOS FL and FL Auto systems as well as the Countess™ II FL Automated Cell Counter. Use the selection guide below to choose the light cube that best matches your dye.

Light cube	Dye	Abs peak	Em peak	Spectral match
GFP	1,8-ANS	372	480	Excellent
	4-methylumbelliferone	360	440	Excellent
	7-amino-4-methylcoumarin (AMC)	351	430	Excellent
	7-hydroxy-4-methylcoumarin	360	449	Excellent
	ABQ	344	445	Excellent
	Acridine	362	462	Excellent
	Alexa Fluor™ 350	346	442	Excellent
	Alexa Fluor 405	401	421	Fair

Light cube	Dye	Abs peak	Em peak	Spectral match
GFP (470/22 Ex; 510/42 Em) Cat. No. AMEP4651	5-Carboxyfluorescein (5-FAM)	492	518	Good
	Acridine orange + DNA	500	526	Good
	Acridine yellow	470	550	Excellent
	ActinGreen™ 488 (Alexa Fluor 488 phalloidin)	496	518	Excellent
	Alexa Fluor 430	432	540	Good
	Alexa Fluor 488	496	518	Excellent
	ATTO 465	453	507	Excellent
	Auramine	460	540	Excellent
	Aurophosphine	470	540	Excellent
	BCECF	499	540	Good
	BODIPY™ 492/515	490	540	Good
	BODIPY 505/515	502	540	Good
	BODIPY FL	502	540	Good
	BTC	464	540	Excellent
	Calcein	494	540	Good
	Calcein Green™ 1	506	540	Fair
		482	540	Excellent

Viability dye	Light cube	Cat. No.
ReadyProbes™ Cell Viability Imaging Kit, Blue/Green	DAPI; GFP	R37609
ReadyProbes™ Cell Viability Imaging Kit, Blue/Red	DAPI; RFP or Texas Red™	R37610
LIVE/DEAD™ Viability/Cytotoxicity Kit	GFP; RFP or Texas Red	L3224
Propidium Iodide ReadyProbes™ Reagent	RFP	R37108
SYTOX™ Green Nucleic Acid Stain	GFP	S7020
SYTOX™ Red Dead Cell Stain	Cy®5	S34859
7-Aminoactinomycin D (7-AAD)	Texas Red or Cy5	A1310

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